# PS300B

Power Supply

## Page Finder

Important Information .....	ii
Waste Electrical and Electronic Equipment (WEEE) .....	vii
Important user information .....	viii
Power supply function and description .....	1
Unpacking .....	3
Specifications.....	4
Parameter overview .....	5
Power supply operation .....	7
Connections .....	8
Operation.....	9
Power Failure Detection (PFD) .....	12
Troubleshooting .....	13
Replacing the fuses .....	14
Care and maintenance .....	15
Technical service and repair .....	15
Ordering information.....	15

## Important Information – English

- If this equipment is used in a manner not specified by Hoefer, Inc. the protection provided by the equipment may be impaired.
- This instrument is designed for indoor laboratory use only.
- Only accessories and parts approved or supplied by Hoefer, Inc. may be used for operating, maintaining, and servicing this product.
- **Warning!** Because this instrument can develop sufficient voltage and current to produce a lethal shock, care must be exercised in its operation.
- This instrument is designed in accordance with the EN61010-1:2001 electrical safety standard. Nevertheless, it should be used only by properly trained operators. Read this entire manual before using the instrument and use only according to the instructions.
- The instrument must always be used with the earth lead of the power cord correctly grounded to earth at the mains outlet.
- Use only undamaged electrical wire and equipment specific for the voltages you will use. All equipment connected to high voltage should be in accordance with EN61010-1:2001.
- Keep the instrument as dry and clean as possible. Wipe regularly with a soft, damp cloth. Let the instrument dry completely before use.
- Do not operate the instrument in extreme humidity (above 80%). Avoid condensation by letting the unit equilibrate to ambient temperature when taking the instrument from a colder to a warmer environment.
- To permit sufficient cooling, ensure that the vents of the instrument are not covered.

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---

## Waste Electrical and Electronic Equipment (WEEE)

English

This symbol indicates that the waste of electrical and electronic equipment must not be disposed as unsorted municipal waste and must be collected separately. Please contact an authorized representative of the manufacturer for information concerning the decommissioning of your equipment.

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---

## Important user information

All users must read this entire manual to fully understand the safe use of the PS300B.

**Please read this entire manual to fully understand the safe and effective use of this product.**

If you have any comments on this manual, please send them to us at:

**Hoefer, Inc.**  
**84 October Hill Road**  
**Holliston, MA 01746**  
**support@hoeferinc.com**

Hoefer, Inc. reserves the right to make changes in the specifications without prior notice.

**WARNING!** The warning symbol highlights instructions that must be followed to avoid personal injury. It is important not to continue until all stated conditions are met and clearly understood. Read the instructions to avoid hazardous conditions.

**CAUTION!** The caution symbol highlights instructions that must be followed to avoid damage to the product or other equipment. It is important not to continue until all stated conditions are met and clearly understood.

**Note:** Note is used to indicate information important for trouble-free and optimal use of the product.

### CE certification

This product meets the requirements of applicable CE-directives. A copy of the corresponding Declaration of Conformity is available on request.

The CE symbol and corresponding declaration of conformity is valid for the instrument when it is:

- used in a laboratory location, and
- used as a stand-alone unit, or
- connected to other CE-marked Hoefer, Inc. instruments, or
- connected to other products recommended or described in this manual, and
- used in the same state as it was delivered from Hoefer, Inc. except for alterations described in this manual.

---

## Power supply function and description

**Note:** The four pairs of output jacks always operate at the same voltage.


The sum of the current of all four outputs is displayed. This value can never exceed 500 mA.

The Hoefer® PS300B power supply can be used with many electrophoresis applications, such as running standard PAGE gels, multiple PAGE minigels, horizontal agarose gels, semidry transfer and minigel tank transfers. The LED display reports power supply status. Parameters are set by pressing UP/DOWN arrow buttons in combination with a SELECT button. Four pairs of output jacks accept 4 mm plugs, which operate in parallel. The output jacks are color coded, recessed for safety, and accept 4 mm shielded (banana) plugs that come standard with most Hoefer equipment.

The user interface consists of a three digit, LED readout, and five buttons on a membrane keypad. Diagrams of the front and rear panel are shown on the following pages.

The PS300B operates in constant voltage or constant current modes. Power output is controlled by setting a maximum value for voltage (up to 300 V), or current (up to 500 mA). The maximum output power at any time is 90 Watts. Run duration can be continuous or programmed to a maximum duration of 999 minutes.


**Fig 1.** Front view of PS300B Power Supply.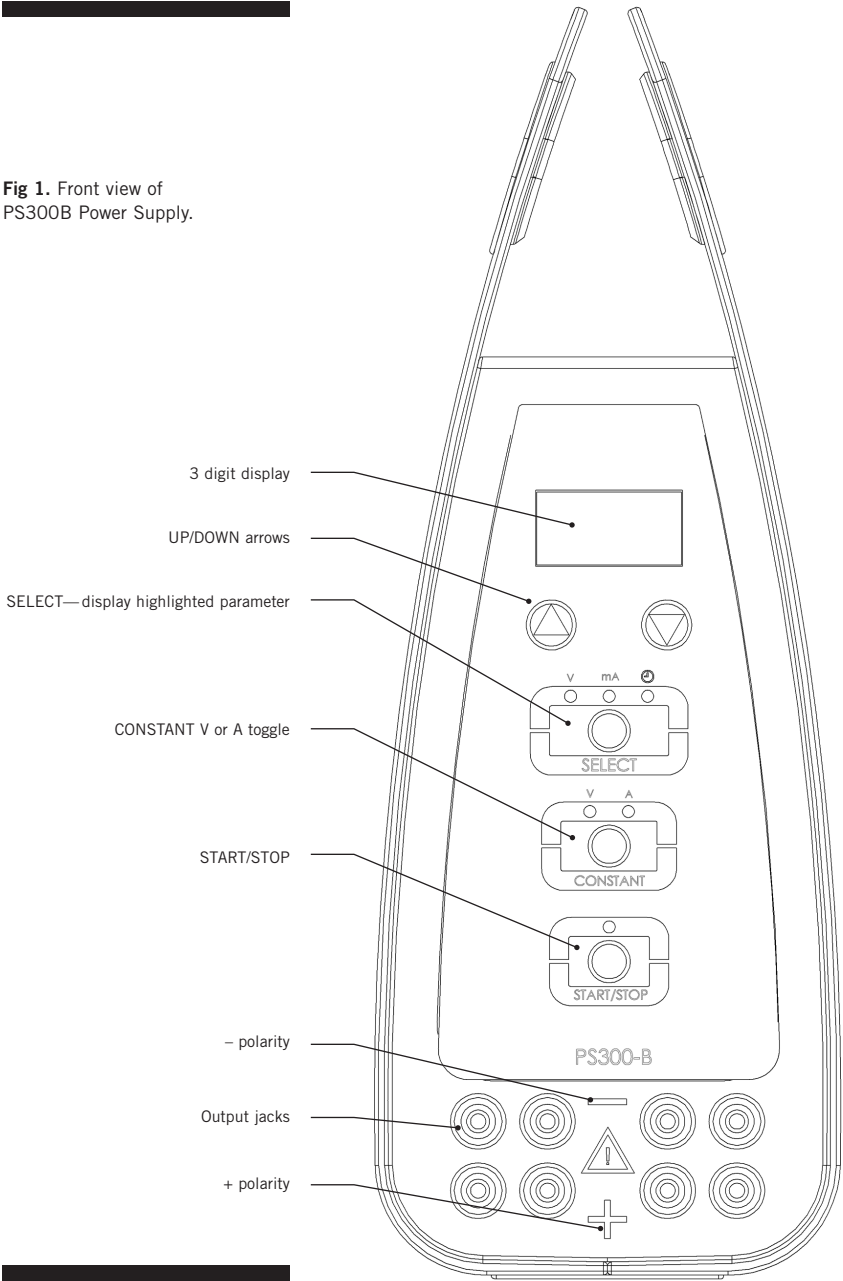

# Unpacking

Unwrap all packages carefully and compare contents with the packing list, making sure all items arrived. If any part is missing, contact your local Hoefer, Inc. sales office. Inspect all components for damage that may have occurred while the unit was in transit. If any part appears damaged, contact the carrier immediately. Be sure to keep all packing material for damage claims or to use should it become necessary to return the unit.

**Fig 2.** Rear view of PS300B Power Supply.

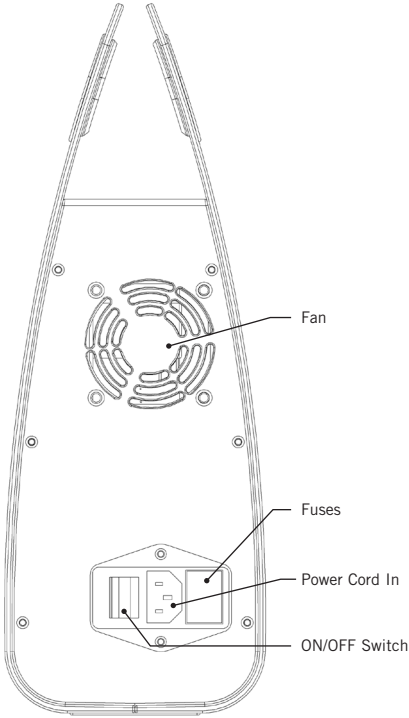## Specifications

<b>Power requirement</b>	Frequency, 50–60 Hz Line voltage, 100–240 VAC 144 VA
<b>User interface</b>	3 digit, red LED membrane keypad
<b>Output</b>	10–300 V, 1 V increment 4–500 mA, 1 mA increment 90 W maximum 4 pairs output jacks in parallel
<b>Output connector</b>	4-mm plugs
<b>Timer</b>	1–999 min, 1 min increment
<b>Environmental operating conditions</b>	Indoor use: 4–40 °C Relative humidity: 0–80% noncondensing Altitude: < 2000 m Pollution degree: 2 Installation category: II
<b>Dimensions (h × w × d)</b>	33.5 × 12.5 × 28 cm 13.2 × 4.9 × 11.0 in
<b>Weight</b>	2.4 kg (5.3 lb)
<b>Safety protection circuits</b>	Short circuit protection Open circuit protection
<b>Product certifications</b>	EN61010-1:2001, EN61326 A3:2003 Class B CE WEEE RoHS

## Parameter overview

### Please note the following power output limitations:

- The maximum power supply output is 90 W, so the maximum current of 500 mA is available at 180 V or less.
- At the maximum voltage setting of 300 V, the current is limited to 300 mA.

**Note:** In most cases, only one parameter can be held constant at a time.

### Power parameters

Because the power output limit is 90 W, and  $W = V \times A$ , the maximum current of 500 mA can be delivered only if voltage is  $\leq 180$  V. Conversely, the maximum voltage of 300 V can be delivered only if the current is  $\leq 300$  mA.

### Programmed limits

Variable	Range	Increment
Voltage	10–300 V	1 volt
Current	4–500 mA	1 milliamp
Time	0–999 minutes	1 minute
Timer = 0 = continuous run		
Timer = 1–999 = timed run in minutes		



### Power supply output limits

The PS300B can be programmed to limit voltage (V) or current (mA). The power supply will operate at the set limits provided the “load,” or power requirement, does not exceed the power supply rating.

“Crossover” occurs when the second parameter limit is reached as system resistance changes (Fig 3). At this point, the second parameter becomes constant, and the previously unchanging parameter adjusts downward as system resistance changes. (System resistance is affected by discontinuous buffers, changes in temperature, and buffer depletion.) When the power supply is delivering power, the light above the CONSTANT button indicates which parameter is being held constant.

Fig 3. Power supply crossover.

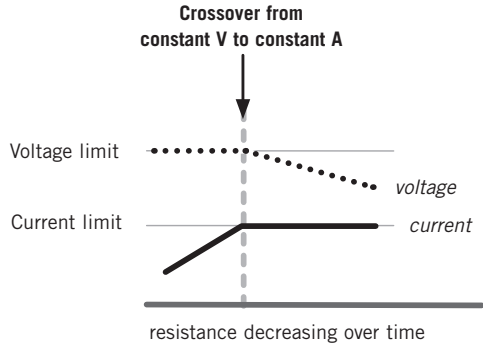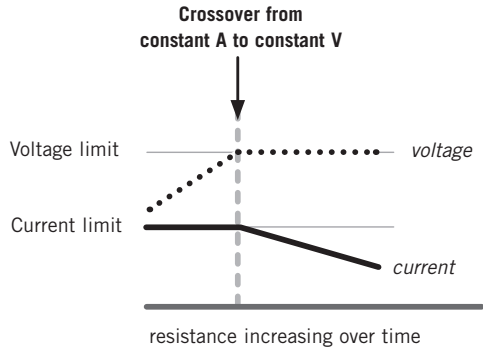There are some conditions when the power supply will operate at the upper limit of 90 W of constant power. When this occurs, both the constant V and the constant mA lights will be lit at the same time.

---

# Power supply operation

**WARNING!** Use only one hand when making or breaking a connection to avoid making a complete circuit across your chest. Also, while connecting or disconnecting the leads, do not allow any part of your body to contact a grounded surface. Make sure your other hand is not touching anything that grounds you.

**WARNING!** If you are in the middle of a run and wish to connect or disconnect the leads, always turn the power supply off (press START/STOP). Wait for the display to read OFF, and the high voltage light to turn off.

## Recommended leads:

- 4 mm (banana) plugs, fixed insulating sheaths, which are found on most Hoefer equipment.

## Will also accept:\*

- 4 mm (banana) plugs, unsheathed or retractable sheaths
- Stackable plugs

\*Exposed metal surfaces on voltage leads can potentially lead to electrical shocks. If using these types of voltage leads, the following safety tips will reduce the risk of electrical shock.

1. Turn off the power supply before connecting or disconnecting the voltage leads.
2. Keep fingers away from metal components when connecting or disconnecting the voltage leads.

# Connections

1

**Connect the mains power cord.** Select the appropriate mains power cord for your region, and connect it to the power cord receptacle on the rear panel and to a suitable grounded AC power outlet. Place the power supply so that the fan on the back panel and the vents on the bottom are not obstructed.


2

**Connect the HV power leads.** Connect the apparatus to the power supply by plugging the lead connectors into one pair of the recessed output jacks (Fig 4).

3

Turn on the main power switch on the rear panel. The front panel should be illuminated.

**Fig 4.** The output jacks are recessed and color coded.

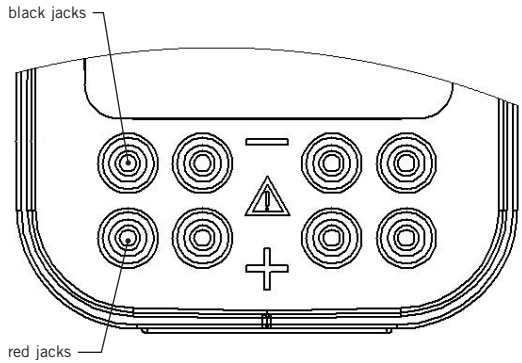# Operation

1

---

**Select constant mode.** Press the CONSTANT button to toggle between constant voltage and constant current mode. The selection is indicated by a green light.

2

---

**Adjust limit.** Press the UP and DOWN buttons to set the constant parameter to the desired value.

3

---

**Check other run parameter.** Press the SELECT button to check the set value of the other (non-constant) run parameter. This parameter is set by default to the maximum output of the power supply. This is the normal setting for many electrophoretic separations.

*Optional:* If there is another upper limit desired for this non-constant parameter, use the UP and DOWN buttons to set it to the desired value. The power supply can then operate in “Crossover” mode as explained on page 5.

4

---

**Set the timer (optional).** Press the SELECT button to move the green indicator light to under the clock symbol.

**Continuous run.** Set the timer at “0” to start a continuous run. The power supply will operate indefinitely, until stopped manually. The timer will start counting up, indicating the length of the run in minutes.

**Timed run.** Press the UP and DOWN buttons to set the timer to the desired run time, from 1 to 999 minutes. The timer counts down during the run, indicating the time remaining.

**Note:** When changing timer between “0” and “1”, a message “PF1” or “PF0” will briefly be displayed, indicating that the Power Failure Detection Feature is either on (PF1) or off (PF0). See page 12 for more information on Power Failure Detection.

**Note:** The power supply slowly increases to the set value when starting a run.

The constant parameter can be adjusted during a run.

## 5

**Start the run.** Press the START/STOP button. The green light above the START/STOP button indicates that voltage and current are being applied to the output jacks. The green light above the CONSTANT button indicates which parameter (V or A) is limiting, and constant. The LED display shows the numerical value of the actual V or A output, or the time (press the SELECT button to toggle between the values).

## 6

**In the case of a continuous run, press START/STOP to terminate the run.** The LED display reads "OFF".

A timed run automatically stops with a warning beep. The word "OFF", and the conditions at the end of the run are flashed on the display so that the user can record the end conditions.

Pushing the SELECT button will return the power supply to the last run settings.

Pressing the CONSTANT button will reset the power supply.

**Example: Set up conditions for running 2 minigels, each at a constant current of 30 mA.**

- 1 Press the CONSTANT button to position the indicator light under "A".
- 2 Press the UP and DOWN buttons to set the LED display value 60. (2 minigels  $\times$  30 mA each).
- 3 Press the SELECT button to position the indicator light under "V". Observe that the volt limit has been automatically set to the maximum value of 300 V.

*Setting the voltage limit to the maximum setting ensures that the entire separation is run at 60 mA for as long as possible. Should the buffer or gel lose too much conductivity during the run, the operational 300 V or 90 W limits may be reached, at which point "crossover" automatically occurs and V becomes the limiting variable. The current would slowly drop below 60 mA at this point.*

- 
- 4 Press the SELECT button to position the indicator light under the Clock symbol. If desired, use the UP button to set the timer to a desired duration. If the run is continuous, set the timer value to “0”.
  - 5 Press the START/STOP button to start the run.

**Example: Set maximum mA (30 mA per gel) and maximum V (180 V) settings, with automatic crossover.**

**Note:** If this crossover condition occurs and is not desired, change the voltage or current setting to the power supply's operational maximum (500 mA or 300 V).

*The power supply can automatically crossover between two different maximum settings, if desired. This will be illustrated below.*

- 1 Press the CONSTANT button to position the indicator light under “A”.
- 2 Press the UP and DOWN buttons to increase the LED display value to 60 (2 minigels × 30 mA each).
- 3 Press the SELECT button to position the indicator light under “V”. Observe that the volt limit has been automatically set to the maximum value of 300 V. Press the UP and DOWN buttons to adjust this value down to 180 V.
- 4 Press the SELECT button to position the indicator light under the Clock symbol. If desired, use the UP button to set the timer to a desired duration. If the run is continuous, set the timer value to “0”.
- 5 Press the START/STOP button to start the run.

*Both Voltage and Current are monitored during the run. The power supply initially outputs 60 mA constant current (2 minigels × 30 mA each). The voltage output will be <180 V, but may rise during the electrophoresis run, owing to gradual depletion of the buffer during the run. If the voltage never exceeds 180 V during the run, then the gels will run at 60 mA for the entire run. If the voltage increases to 180 V, then the power supply will automatically “crossover” to constant voltage mode, in order to keep voltage from exceeding 180 V. Now, output current will slowly decrease below the set limit of 60 mA because the 180 V set limit is now the limiting value in the circuit.*

## Power Failure Detection (PFD)

The Power Failure Detection feature will save the time, voltage and current settings in the event of a power outage, and automatically continue the run when the power is restored.

To alert the user that a power failure occurred during the run, the error message “E6” flashes at the end of the run, instead of the usual “OFF” message. The “E6” error message means that despite the power outage, the power supply successfully supplied the programmed voltage and current for the entire length of time that was programmed. The time shown at the end of the run is the time the gel ran since power was restored, not the total time of the run.

**The default settings for Power Failure Detection are:**

Run	Time Setting	PFD	Code
Timed Run	1–999 minutes	ON	PF1
Continuous Run	0	OFF	PFO

**The Power Failure Detection feature can be deactivated during a timed run.**

- 1 Press the SELECT button to move the indicator light under the Clock symbol.
- 2 Press the UP and DOWN buttons to enter in the run time in minutes.
- 3 Press both the UP and DOWN buttons at the same time. The display will briefly show “PFO”, indicating that the PFD mode is disabled.
- 4 Press both the UP and DOWN buttons at the same time again. The display will show “PF1”, indicating that the PFD mode is enabled.

# Troubleshooting

The power supply has built-in error messages to help troubleshoot problems with the electrical circuit. The error messages, and possible corrective actions, are listed below:

Error code	Explanation	Possible solutions
E1	<b>No Load Detected.</b> The voltage leads to the electrophoresis cell are not attached. The current load is below 4 mA, the lower limit of the power supply.	Check all connections. Verify that buffer volumes are sufficient to contact all electrode wires and that buffer is contacting both surfaces of the gel. A constant volt run with low volt setting may result in a current below the 4 mA limit. Press any button* to clear the error code.
E3	<b>Change in Load Resistance.</b> Open circuit was detected at some point after a successful start. Buffer is leaking from the cells. There is a loose electrical connection in the circuit.	Check and correct all potential resistance problems. Check buffer volumes. Check all voltage leads. Check the fit of power supply adapters, if used. The output current may have fallen below the 4 mA limit of the power supply. Press any button* to clear the error code.
E4	<b>Power Failure</b> during a continuous run (no timer being used)	Restart the run, if desired. Press any button* to clear the error code.
E5	<b>Power Failure</b> during a timed run with PFD option disabled. Run was not completed.	Restart the run, if desired. Press any button* to clear the error code. See page 12 for information on the PFD option.
E6	<b>Power Failure</b> during a timed run with PFD option enabled. Despite the power outage the run was completed successfully.	Press any button* to clear the error code. See page 12 for information on the PFD option.
E7	<b>Short Circuit</b> Load current exceeded 500 mA. The voltage load is below 10 V, the lower limit of the power supply.	Check for any short circuit or evidence of electrical arcing. Buffer concentration too high. A constant current run with low current may result in a voltage below the 10 V limit.

\*Press any button, except the START/STOP button, to clear the Error Message.

## Other Problems:


**No display and/or fan not operating:** Check that the power supply is plugged into a working receptacle. Check that the main power switch is on. Check fuses.

**CAUTION!** Turn the mains power supply switch off and detach the power cord before replacing input mains fuses!

## Replacing the fuses

- Unplug the power cord from the power supply.
- Locate the fuse compartment on the rear panel.
- Use a screwdriver to pry out the fuse compartment. There are two fuses in the compartment. Check and replace any blown fuses with a fuse of the same size and ratings.
- FUSES (2); 250 V 3.15 A 5 × 20 mm fast blow.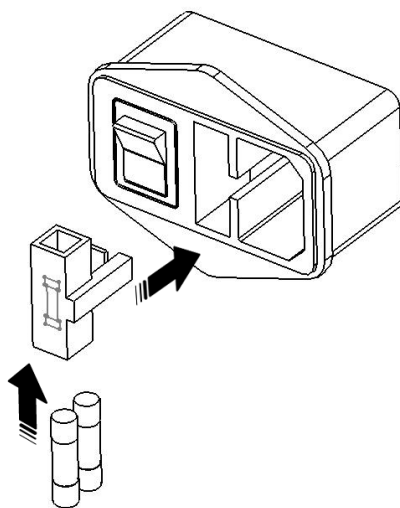

## Care and maintenance

- Turn the mains power switch off and unplug the power cord before cleaning.
- Use a soft cloth dampened with water or a mild cleaning solution to clean the cabinet and display.
- If spilled liquids contact the circuit boards, unplug the power supply and allow to dry completely. Call Hoefler, Inc. for advice before using.

## Technical service and repair

Hoefler, Inc. offers complete technical support for all of our products. If you have any questions about how to use this product, or would like to arrange to repair it, please call or fax your local Hoefler, Inc. representative.

## Ordering information

	Qty.	Code no.
PS300B Power Supply, universal voltage	1	PS300B
Power Cord, US	1	PSCORD-115V
Power Cord, Euro	1	PSCORD-230V
User Manual	1	PS300B-IM
Fuse 250 V 3.15 A FB 5 × 20 mm	pack/5	PSF3A-FB-5X20

**Hoefer, Inc.**

84 October Hill Road  
Holliston, MA 01746

Toll Free: 1-800-227-4750

Phone: 1-508-893-8999

Fax: 1-508-893-0176

E-mail: [support@hoeferinc.com](mailto:support@hoeferinc.com)

Web: [www.hoeferinc.com](http://www.hoeferinc.com)

Hoefer is a registered trademark  
of Hoefer, Inc.

© 2012 Hoefer, Inc. —

All rights reserved.

Printed in the USA.